Delphi-X Observer

Superior imaging






State-of-the-art microscopes for advanced Life Science applications

The ergonomic design and outstanding state-of-the-art objectives make the Delphi-X Observer the ideal microscope for advanced Life Science applications such as cytology and anatomopathology

The 25 mm field of view of the eyepieces and the plan apochromatic objectives enable observations with perfect color rendering at high resolving power

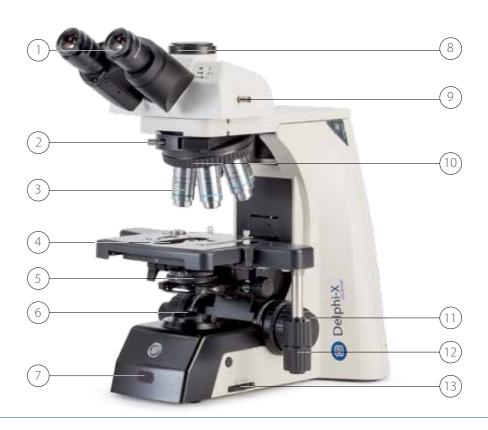

FEATURES

- 1 | DIN Super Wide field SWF 10x/25 mm, Ø 30 mm
- 2 | For slider with 1 lambda retardation plate for observation of monosodium crystal deposition (gout) and for DIC slot (DIC ready)
- **3** Infinity corrected 60 mm achromatic Plan, Plan
 Phase or Plan Semi-Apochromatic EIS objectives
- 4 | 300/190 x 152 mm stage with 78 x 32 mm integrated right-handed mechanical stage with scratch-resistant Gorilla glass
- **5** | Height adjustable N.A. 0.90 /1.25 Abbe swing-out condenser

- 6 | 3 W LED or 100 W halogen light source
- 7 | iCare sensor for energy saving
- 8 | Universal photo tube
- **9** | 3-Position beamsplitter (100:0 / 80:20 / 0:100)
- **10** | Revolving sextuple reversed nosepiece on ball-bearings
- **11** | Precise coarse and fine adjustments
- **12** | Low stage adjustments
- **13** | Full Köhler diaphragm

OPTIONAL

- EWF 10x/22 mm eyepieces, Ø 30 mm
- Left-handed stage
- Stage with scratch-resistant Sapphire glass
- Ergonomic tilting head
- Face-to-face dual head
- Multi head with 2, 3,5 or 10 heads

Welcome to the new world of advanced microscopy

With 25 mm field of view, unsurpassed 60 mm EIS objectives, sextuple revolver and a large rackless stage with scratch-resistant Gorilla glass, the Delphi-X Observer offers the absolute best available Life Science microscope on the market. Long working sessions will not be a problem due to the ergonomic design, low stage controls for either left or right-handed users and an optional ergonomic tilting head to match each user requirement



EYEPIECES

- DIN Super Wide field SWF 10x/25 mm, Ø 30 mm eyepieces
- DIN Super Wide Field EWF 10x/22 mm, Ø 30 mm eyepieces also available on special request

HEAD

- Siedentopf trinocular with 30° inclined tubes
- Interpupillary distance adjustable between 47 and 78 mm
- 3-Position beamsplitter (100:0 / 80:20 / 0:100)
- Diopter adjustment on both eyepieces
- · Optional: ergonomic tilting head
- Optional: face-to-face dual head
- Optional: multi head system with 2, 3, 5 or 10 heads

NOSEPIECE

· Revolving sextuple reversed nosepiece on ball-bearings



Optional: ergonomic tilting head



For slider with 1λ retardation plate for observation of monosodium crystal deposition (gout) and for DIC slot (DIC ready)



The sextuple revolver allows users to choose six 60 mm EIS objectives with the highest possible resolving power and maximum contrast Designed to be superior in accuracy and color reproduction, the EIS objectives are virtually free of aberrations 6 | @euromex®

Observe details you have never seen before

The new design of infinity corrected Plan, Plan phase and Semi-Apochromatic 60 mm EIS objectives produce images with the highest possible resolving power and contrast. Virtually free of image distortion the image quality is absolutely superb and will let you observe details you have never seen before.

With a field of view up to 25 mm, the Delphi-X Observer is suitable for the most demanding applications



DX.1153-PLi

ENHANCED INFINITY SYSTEM (EIS)

With 25 mm field of view, unsurpassed 60 mm Enhanced Infinity System (EIS) objectives, sextuple revolver and a large rackless stage with scratch-resistant Gorilla glass the Delphi-X Observer offers the absolute best available Life Science microscope on the market. Long working sessions will not be a problem due to the ergonomic design, low stage controls for either left or right handed users and an optional ergonomic tilting head to match each user requirement

All optics are anti-fungus treated and anti-reflection coated for maximum light throughput. The S40x, S50x, S60x and S100x objectives are spring loaded

OBJECTIVES

(ANATOMO- AND HISTOPATHOLOGY LABORATORIES)

 Infinity corrected Plan Semi-Apochromatic EIS 4x/0.13 WD 16.5 mm, 10x/0.30 WD 8.1 mm, 20x/0.50 WD 2.1 mm, S40x/0.75 WD 0.70 mm, S100x/1.30 (oil immersion) 0.15 mm objectives

OBJECTIVES

(CLINICAL ROUTINE LABORATORIES)

- Infinity corrected Plan Achromatic EIS 4x/0.10 WD 30 mm, 10x/0.25 WD 10.2 mm, 20x/0.40 WD 12.0 mm, S40x/0.65 WD 0.7 mm, S100x/1.25 (oil immersion) WD 0.2 mm objectives
- Optional 2x/0.06 WD 7.5 mm, S50x/0.95 (oil immersion) WD 0.19 mm and S60x/0.80 WD 0.30 mm

OBJECTIVES

(PHASE CONTRAST)

 Infinity corrected Plan Phase EIS 10x/0.25 WD 10.2 mm, 20x/0.40 WD 12 mm, S40x/0.65 WD 0.7 mm, S100x/1.25 (oil immersion). WD 0.2 mm objectives

WD = Working distance

OBJECTIVES

Plan PLi EIS	2x/0.06 ⁽¹⁾	4x/0.10	10x/0.25	20x/0.40	S40x/0.65	S50x/0.95 oil (1)	S60x/0.80 (1)	S100x/1.25 oil	S100x/1.25 oil iris (1)
Plan Phase PLPHi EIS			10x/0.26	20x/0.41	S40x/0.66			S100x/1.25 oil	
Plan Semi-Apochromatic		4x/0.13	10x/0.30	20x/0.50	S40x/0.75			S100x/1.30 oil	

(1) Optional

Designed with ergonomics in mind

The Delphi-X Observer's design provides more comfort and convenience for professional microscopists. The double fine and coarse focus control knobs can be switched from left to right as per user preference. The anti-slip coated adjustment knobs provide maximum user comfort during displacement of the sample. Switching of the neutral density filters, setting the Köhler diaphragm and adjusting light intensity can all be controlled without lifting your hands to prevent fatigue



Optional: mechanical stage with scratch-resistant Sapphire glass

STAGE

- 300/190 x 152 mm stage with 78 x 32 mm integrated right-handed mechanical stage
- · Left-handed mechanical stage available on request
- Standard supplied with robust and scratch-resistant Gorilla glass
- · Optional Sapphire glass available on request

FOCUSING

- Coaxial coarse and fine adjustment, 100 graduations, 1 μm precision, 100 µm per rotation, total travel approximately 35 mm
- Supplied with an adjustable rack stop to prevent damage to sample and objectives
- ${\boldsymbol{\cdot}}$ The coarse adjustments are equipped with friction control

CONDENSER FOR BRIGHTFIELD

• In height adjustable N.A. 0.90/1.25 Abbe swing-out condenser

CONDENSER FOR PHASE CONTRAST

• In height adjustable N.A. 1.25 Zernike phase contrast condenser with phase annuli for 10/20/S40x and S100x-oil objectives







The carrying grip at the back of the microscope ensures safe transportation of the microscope and the integrated tool & holder make sure the right tool is always available

DX.1153-PLi

DIASCOPIC ILLUMINATIONS

Two illumination versions are available for the Delphi-X Observer microscope:

- Diascopic intensity adjustable 3 W NeoLED $\!^{\mathrm{m}}$ illumination with internal 100-240 Vac power supply. The diascopic NeoLED™ illumination comes with 1 neutral density filter
- Diascopic intensity adjustable 100 W halogen illumination with internal 100-240 Vac power supply. The diascopic halogen illumination comes with 2 push-in/push-out neutral density filters for smooth attenuation of the light intensity for all kind of samples



iCare sensor for energy saving



Filter selection button

NEOLED

The innovative NeoLED $^{\mathrm{IM}}$ design is a combination of a custom LED and a specially designed thin lens with a short focal length in order to obtain three main benefits:

- More oblique light from the LED light source can be captured, which increases the light output significantly
- Less energy is required to achieve this level of light intensity
- The larger aperture of NeoLED™ allows the optical system of the microscope to produce images at higher resolutions, very close to the theoretical diffraction limit of the optics

ICARE SENSOR

The unique iCare Sensor is developed to avoid unnecessary loss of energy. The illumination of the microscope automatically switches off shortly after microscopists step away from their position

PACKAGING

Supplied with power cord, dust cover, spare fuse, user manual and a bottle of 5 ml immersion oil

Simultaneous observation of specimens

Hospitals, medical centers and universities highly appreciate the unique Delphi-X Observer multi head systems. The high quality image can simultaneously be observed up to 10 colleagues or students at the same time

All systems are equipped with a dual color (green or red) laser pointer that allows the main user to point out areas of interest across the entire field of views of all heads by using the built-in joystick



MULTI HEAD SYSTEM

The Delphi-X Observer can be extended with a unique multi head system for in total 2, 3, 5 or 10 heads allowing simultaneous observation of specimens

The central attachment is equipped with a trinocular head and with a joystick to move a green or red laser pointer across the entire field of views of all heads. Configurations with one, two, four or nine additional binocular heads, 30° inclined tubes and infinity corrected IOS objectives.

The trinocular head is supplied with extended wide field EWF 10x/25 mm eyepieces. Each extra binocular head is supplied with extended wide field plan EWF 10x/25 mm eyepieces except for the model with 9 binocular heads (10 users) that is supplied with EWF 10x/22 mm eyepieces

FACE-TO-FACE DUAL HEAD SYSTEM

The Delphi-X Observer can also be equipped with a face-to-face trinocular head for 2 users allowing simultaneous observation

Supplied with 30° inclined tubes and plan infinity corrected IOS objectives. The central attachment is equipped with a joystick to move a green or red laser

pointer across the entire field of views of the 2 heads. The trinocular head is supplied with extended wide field EWF 10x/25 mm eyepieces and the second binocular head is supplied with extended wide field EWF 10x/22 mm

Delphi-X Observer face-to-face dual head system





MODELS WITH TRINOCULAR HEAD

	Trinocular	SWF 10x/25 mm eyepieces	Sextuple nosepiece	Plan 4/10/20/S40/S100 oil EIS objectives	Plan Semi-Apochromatic 4/10/20/S40/S100 oil EIS objectives	Plan Phase 10/20/S40/S100 oil EIS objectives	3 W NeoLED Köhler	100 W 24 V halogen Köhler
DX.1153-PLi	•	•	•	•			•	
DX.1153-PLFi	•	•	•		•			
DX.1153-PLPHi	•	•	•			•	•	
DX.2153-PLi	•	•	•	•				•
DX.2153-PLFi		•						
DX 2153-PI PHi								

MODELS WITH ERGONOMIC TRINOCULAR TILTING HEAD

	Trinocular	SWF 10x/25 mm eyepieces	Sextuple nosepiece	Plan 4/10/20/S40/S100 oil EIS objectives	Plan Semi-Apochromatic 4/10/20/S40/S100 oil EIS objectives	Plan Phase 10/20/S40/S100 oil EIS objectives	3 W NeoLED Köhler	100 W 24 V halogen Köhler
DX.1158-PLi	•	•	•	•			•	
DX.1158-PLFi	•	•	•		•		•	
DX.1158-PLPHi	•	•	•			•	•	
DX.2158-PLi	•	•	•	•				
DX.2158-PLFi	•	•	•					•
DX.2158-PLPHi	•	•	•					

MULTI HEAD SYSTEMS

	Total users	Trinocular SWF 10x/25 mm eyepieces	Binocular SWF 10x/25 mm eyepieces	Binocular SWF 10x/22 mm eyepieces	Plan 4/10/20/S40/S100 oil EIS objectives	Plan Semi-Apochromatic 4/10/20/S40/S100 oil EIS objectives	Plan Phase 10/20/S40/S100 oil EIS objectives	3 W NeoLED Köhler	10 W NeoLED Köhler	100 W 24 V halogen Köhler	iCare sensor
DX.1156-PLi/2	2	1	1	-	•			•			•
DX.1156-PLFi/2	2	1	1	-		•		•			•
DX.1156-PLPHi/2	2	1	1	-			•	•			•
DX.2156-PLi/2	2	1	1	-	•					•	•
DX.2156-PLFi/2	2	1	1	-		•				•	•
DX.2156-PLPHi/2	2	1	1	-			•			•	•
DX.1156-PLi/3	3	1	2	-	•			•			•
DX.1156-PLFi/3	3	1	2	-		•		•			•
DX.1156-PLPHi/3	3	1	2	-			•	•			•
DX.2156-PLi/3	3	1	2	-	•						•
DX.2156-PLFi/3	3	1	2	-						•	•
DX.2156-PLPHi/3	3	1	2	-			•			•	•
DX.1156-PLi/5	5	1	4	-	•			•			•
DX.1156-PLFi/5	5	1	4	-		•		•			•
DX.1156-PLPHi/5	5	1	4	-			•	•			•
DX.2156-PLi/5	5	1	4	-	•					•	•
DX.2156-PLFi/5	5	1	4	-		•				•	•
DX.2156-PLPHi/5	5	1	4	-			•			•	•
DX.1156-PLi/10	10	1	-	9	•				•		
DX.1156-PLFi/10	10	1	-	9		•			•		
DX.1156-PLPHi/10	10	1	-	9			•		•		
DX.2156-PLi/10	10	1	-	9	•					•	•
DX.2156-PLFi/10	10	1	-	9		•				•	•
DX.2156-PLPHi/10	10	1	-	9			•			•	•

FACE-TO-FACE DUAL HEAD SYSTEMS

	Total users	Trinocular SWF 10x/25 mm eyepieces	Binocular EWF 10x/22 mm eyepiece	Sextuple nosepiece	Plan 4/10/20/S40/S100 oil EIS objectives	Plan Semi-Apochromatic 4/10/20/S40/S100 oil EIS objectives	Plan Phase 10/20/S40/S100 oil EIS objectives	3 W NeoLED Köhler	100 W 24 V halogen Köhler
DX.1154-PLi	2	1	1		•			•	
DX.1154-PLFi	2	1	1	•		•		•	
DX.1154-PLPHi	2	1	1	•				•	
DX.2154-PLi	2	1	1	•	•				•
DX.2154-PLFi	2	1	1	•		•			•
DX 2154-PI PHi	2	1	1						

Left-handed mechanical stage available on request

14 | @ euromex°

Accessories and spare parts

ALL MO	DELS		
DX.6010	Super wide field SWF 10x/25 mm eyepiece, Ø 30 mm tube	DX.7710	Plan Phase PLPHi 10x/0.25 infinity corrected objective.
DX.6010-C	Super wide field SWF 10x/25 mm eyepiece		Working distance 10.2 mm
	with cross-hair, Ø 30 mm tube	DX.7720	Plan Phase PLPHi 20x/0.40 infinity corrected objective.
DX.6010-N	1 Super wide field SWF 10x/25 mm eyepiece		Working distance 12 mm
	with micrometer, Ø 30 mm tube	DX.7740	Plan Phase PLPHi S40x/0.65 infinity
DX.6010-C	M Super wide field SWF 10x/25 mm eyepiece		corrected objective. Working distance 0.7 mm
	with micrometer and cross-hair, Ø 30 mm tube	DX.7700	Plan Phase PLPHi S100x/1.25 oil immersion
DX.6010-S	G Super wide field SWF 10x/25 mm eyepiece		infinity corrected objective. Working distance 0.2 mm
	with 20 x 20 square grid reticle, Ø 30 mm tube		
		DX.9105	Abbe condenser 0.9/1.25 with swing-out front lens
DX.6210	Super wide field SWF 10x/22 mm eyepiece, Ø 30 mm tube	DX.9126	Zernike phase contrast set
DX.6210-C	Super wide field SWF 10x/22 mm eyepiece		with Plan Phase PLPHi 10/20/S40/S100x objectives,
	with cross-hair, Ø 30 mm tube		phase contrast condenser, telescope and green filter
DX.6210-N	1 Super wide field SWF 10x/22 mm eyepiece	DX.9140	Zernike Phase contrast Abbe condenser with 4 phase annuli
	with micrometer, Ø 30 mm tube	DX.9148	Telescope for phase contrast alignment, Ø 30 mm tube
DX.6210-C	M Super wide field SWF 10x/22 mm eyepiece		
	with micrometer and cross-hair, Ø 30 mm tube	DX.9660	Simple polarization kit,
DX.6210-S	G Super wide field SWF 10x/22 mm eyepiece		consisting of polarizer in slider and analyzer on lamp house
	with 20 x 20 square grid reticle, Ø 30 mm tube	DX.9662	Simple polarization kit
			with 1λ retardation plate for identification of gout crystals
DX.6012	Wide field WF 12.5x/17.5 mm eyepiece, Ø 30 mm tube		(monosodium urate)
DX.6015	Wide field WF 15x/16 mm eyepiece, Ø 30 mm tube		
DX.6020	Wide field WF 20x/12 mm eyepiece, Ø 30 mm tube	DX.9501-I	L G Option for left-handed stage with Gorilla glass.
DX.6099-L	Eyecups for eyepieces		Only with new Delphi-X Observer microscopes (DX.xx5x-xxx/LG
		DX.9504-I	RS Option for right-handed stage with Sapphire glass.
DX.7202	Plan achromatic PLi 2x/0.06 infinity corrected objective.		Only with new Delphi-X Observer microscopes (DX.xx5x-xxx/ RS
	Working distance 7.5 mm	DX.9504-I	LS Option for left-handed stage with Sapphire glass.
DX.7204	Plan achromatic PLi 4x/0.10 infinity corrected objective.		Only with new Delphi-X Observer microscopes (DX.xx5x-xxx/ LS
	Working distance 30 mm		
DX.7210	Plan achromatic PLi 10x/0.25 infinity corrected objective.	DX.9700	Blue filter, Ø 45 mm, fits on lamphouse
	Working distance 10.2 mm	DX.9702	Green filter, Ø 45 mm, fits on lamphouse
DX.7220	Plan achromatic PLi 20x/0.40 infinity corrected objective.	DX.9704	Yellow filter, Ø 45 mm, fits on lamphouse
	Working distance 12 mm		
DX.7240	Plan achromatic PLi S40x/0.65 infinity corrected objective.	DX.9930	NeoLED™ replacement unit
	Working distance 0.7 mm	SL.3680	100 W 24 Vac halogen replacement bulb
DX.7250	Plan achromatic PLi S50x/0.95 oil immersion infinity	SL.1385	100 W HBO mercury vapor lamp
	corrected objective. Working distance 0.19 mm		
DX.7260	Plan achromatic PLi S60x/0.80 infinity corrected objective.	AE.5214	Fuses 500 mA 250 V, 10 pieces
	Working distance 0.3 mm		(For Delphi-X Observer models with LED)
DX.7200	Plan achromatic PLi S100x/1.25 oil immersion	AE.5216	Fuses 5 A 250 V, 10 pieces
	infinity corrected objective. Working distance 0.2 mm		(For Delphi-X Observer models with 100 W halogen)
DX.7200-I	Plan achromatic PLi S100x/1.25 oil immersion		
	infinity corrected objective with built-in iris diaphragm.	AE.5130	Universal Ø 23.2 mm tube adapter with built-in 2x lens
	For darkfield applications. Working distance 0.2 mm		for SLR photo camera with APS-C sensor. Needs T2 adapter
		AE.5025	T2 ring for Nikon D SLR digital camera
DX.7304	Plan Semi-Apochromatic PLi-Apo 4x/0.13	AE.5040	T2 ring for Canon EOS SLR digital camera
	infinity corrected objective. Working distance 16.5 mm		
DX.7310	Plan Semi-Apochromatic PLi-Apo 10x/0.30	PB.5155	Microscope slides 76 x 26 mm, ground edges, 50 pieces
	infinity corrected objective. Working distance 8.1 mm	PB.5165	Cover glasses 18 x 18 mm, thickness 0.13-0.17 mm, 100 pieces
DX.7320	Plan Semi-Apochromatic PLi-Apo 20x/0.50	PB.5168	Cover glasses 22 x 22 mm, thickness 0.13-0.17 mm, 100 pieces
	infinity corrected objective. Working distance 2.1 mm	PB.5245	Lens cleaning paper, 100 sheets per pack
DX.7340	Plan Semi-Apochromatic PLi-Apo S40x/0.75	PB.5255	Immersion oil, n = 1.482 (25 ml)
	infinity corrected objective. Working distance 0.7 mm	PB.5274	Isopropyl alcohol 99% (200 ml)
DV 7200	Di . C A I	PR 5275	Cleaning kit-lens cleaning fluid lint free lens tissue

PB.5275 Cleaning kit: lens cleaning fluid, lint free lens tissue,

brush, air blower, cotton swabs

DX.7300 Plan Semi-Apochromatic PLi-Apo S100x/1.3 oil immersion

infinity corrected objective. Working distance 0.15 mm

Delphi-X Observer Superb in fluorescence

For biological, medical or biotechnology applications, fluorescence microscopes are used to observe living cells or tissues with help of fluorophores, which act as biomarkers in the cells or tissues to be observed. When exposed to a fluorescence light source, excitation energy is absorbed and starts emitting so-called emission light that can be observed through the eyepieces of the microscope




The state-of-the-art optics and mechanical components make the Delphi-X Observer suitable for routine and advanced fluorescence applications in Life Science, Biotechnical and Medical fields

ILLUMINATION FOR FLUORESCENCE

The Delphi-X Observer for fluorescence is equipped with a 6-position rotating attachment with 100 W mercury vapor light source for broad powerful light spectrum, UV protection shield, integrated safety shutter and a 100-240 Vac external power supply with indication of the lifetime of the 100 W HBO lamp

The 6-position rotating attachment can be supplied with general purpose fluorescence filters for routine applications for blue, green, violet and ultra-violet excitation or with advanced fluorescence filters optimized for specific fluorophores All Delphi-X Observer fluorescence filter sets are supplied with state-of-the-art sputtered and hard coated band-pass excitation filters, high-efficiency dichroic mirrors with very high slopes for superior spectral separation and high blocking long-pass emission filters, all mounted

in anodized metal cubes

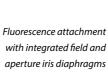







Filter set for blue excitation mounted in filter cube for 6-position fluorescence attachment

DX.9745 / DX.9749

SELECTION OF **FLUORESCENCE FILTER SETS**

Choosing the appropriate filter set is a very important step before ordering your filters. It is essential to match the spectral characteristics of excitation and emission filters and dichroic mirror with the fluorophores used in your fluorescence microscopy application

The most important characteristics for the selection of fluorescence filter sets are:

- The separation between excitation and emission wavelengths
- The slope of the dichroic mirror
- The optical density of the emission filters
- The type of emission filter: longpass or bandpass in terms of color discrimination
- The transmission percentage of the excitation and emission filters
- The type of coating applied to the filters in terms of durability

STANDARD FLUORESCENCE FILTER SETS

Our standard fluorescence filter sets consists out of standard sputtered bandpass filter sets, high sharp slope dichroic mirrors for excellent separation and emission filters with a high transmission level above 90%. These filter sets are offered at an affordable price and have a high hardness and excellent durability

These filter sets are perfect for routine fluorescence and for most applications requiring precise color discrimination and excellent out-of-band blocking. For applications requiring the highest possible signal intensities, optimized and optimized premium filter sets are recommended

OPTIMIZED AND OPTIMIZED PREMIUM FLUORESCENCE FILTER SETS

These filters sets feature the highest throughput sputter coatings with maximum transmission levels. The filters have very steep transitions and very high optical density suitable for modern fluorophores with short Stokes shifts (difference between absorption and emission maxima) and very little spectral separation between excitation and emission bands

The high hardness and excellent durability of the coatings enables these fluorescence filter sets to be used in high heat and humidity

6-POSITION FLUORESCENCE ROTATING ATTACHMENT

The 6-position rotating attachment can be supplied with standard fluorescence filters for routine applications for blue, green, violet and ultra-violet excitation or advanced fluorescence filters optimized for specific fluorophore(s)

FLUORESCENCE MODELS

	Trinocular	SWF 10x/25 mm eyepieces	Diopter adjustable (both sides)	Plan 4/10/20/40/ S100x EIS objectives	Plan Fluarex 4/10/20/40/ S100x objectives	100 W HBO fluorescence	Filter sets (blue&green)	3 W NeoLED	100 W 24 V halogen
DX.3153-PLi	•	•	•	•		•	•	•	
DX.3153-PLFi	•	•	•		•	•	•	•	
DX.3153-PLi/HAL	•	•	•	•		•	•		•
DV 3153 DI EI/HAI	_	_			_	_	_		_

Left-handed mechanical stage available on request

Accessories and spare parts

STANDARD FLUORESCENCE FILTER SETS

Needs empty filter block

DX.9745 Filter set for blue excitation,

FX: 460-490 nm BP. DM: 500 nm. FM: 520 nm LP

DX.9746 Filter set for green excitation,

EX: 510-550 nm BP, DM: 570 nm, EM: 590 nm LP

DX.9747 Filter set for violet excitation.

FX: 400-410 nm RP. DM: 455 nm. FM: 460 nm LP DX.9748 Filter set for ultra-violet excitation,

EX: 330-385 nm BP, DM: 400 nm, EM: 420 nm LP

OPTIMIZED FLUORESCENCE FILTER SETS

Needs empty filter block

DX.9750.39001

Filter set EX435 (BP20) DM455 (LP) EM480 (BP30) AmCyan, Atto 425, Cerulean, CFP, ECFP fluorophore

DX.9750.39002

Filter set EX480 (BP30) DM505(LP) EM535 (BP40) Acridine Orange + DNA, Alexa Fluor 488™, Atto 488, Azami Green, BODIPY FL/pH7.2, Calcein, Calcium Green[™]-1, Cy2[™], DiO, DyLight 488, EGFP, Emerald GFP, FAM, FITC, Fluo-4, GFP, LysoTracker Green/pH 5.2, MitoTracker Green, FM/MeOH, mWasabi, Oregon, Green™ 488, SYBR® Green I, SYTO 9/DNA, ZsGreen fluorophore

DX.9750.39003

Filter set EX495 (BP20) DM515 (LP) EM540 (BP30) Alexa Fluor 514[™], Citrine, EYFP/pH 7, F/IAsH-CCPFCC, mCitrine, Oregon Green™ 514, Rhodamine 123, Topaz

DX.9750.39004

Filter set EX540 (BP25) DM565 (LP) EM605 (BP55) Alexa Fluor 546[™], Alexa Fluor 555[™], Atto 550, CAL Fluor® Red 590, Cy3[™], Dil, DsRed, DyLight 549, MitoTracker Orange/MeOH, Resorufin, Rhod-2, TagRFP, TAMRA, tdTomato Tetramethylrhodamine isothiocyanate, TRITC

DX.9750.39005

Filter set EX540 (BP25) DM565 (LP) EM 620 (BP60) TRITC-red-shifted, Ethidium homidimer-1/DNA, Propidium Iodide

DX.9750.39007

Filter set EX620 (BP50) DM655 (LP) EM 690 (BP50) Alexa Fluor 647™, Allophycocyanin (APC), Atto 647N, Cy5™, DiD, Drag5, DayLight 649, MitoTracker Deep Red 633/MeOH, Nile Blue, Quasar® 670, SYTO® 60, TO-PRO™-3

DX.9750.39008

Filter set EX420 (BP40) DM455 (LP) EM 535 (BP40) Lucifer Yellow/ Tetracycline/AlexaFluor 430

DX.9750.39009

Filter set EX480 (BP30) DM505 (LP) EM 620 (BP60) for Acridine Orange + RNA, D i-8-ANEPP S non-ratiometric, DiA, Ethidium, Bromide, FM™ 1-43, LysoTracker Yellow HCK-123, Pulsar™ 650

OPTIMIZED PREMIUM FLUORESCENCE FILTER SETS

Needs empty filter block

DX.9750.0401

Premium filter set EX440 (BP21), DM455 (LP), EM480 (BP30) CFP, Cerulean, CyPet, SpectrumAqua

DX.9750.0403

Premium filter set EX380 (BP50), DM410 (LP), EM450 (BP60) BFP, BFP2, DAPI, Alexa Fluor 350, Hoechst 33342 & 33258, BFP

DX.9750.0404

Premium filter set EX470 (BP40), DM500 (LP), EM535 (BP50) Alexa Fluor 488, CoralHue Azami Green, Cy2, eGFP, FITC, GFP (sapphire), Emerald

DX.9750.0407

Premium filter set EX635 (BP30), DM660 (LP), EM710 (BP80) Cy5, Alexa Fluor647. Also compatible with TO-PRO-3, TOTO-3, Kodak, X-Sight 650

DX.9750.0412

Premium filter set EX500 (BP25), DM252 (LP), EM545 (BP35) Acridine Orange (+DNA), Rhodamine Green, CL NERF, (hi pH), YFP. Also compatible with Alexa Fluor 532

Premium filter set EX560 (BP55), DM505 (LP), EM645 (BP75) Texas Red, Texas Red-X, CryptoLight CF-6. Also compatible with Cy3.5, HcRed, mRaspberry, MitoTracker Red

OTHER

DX.9200-6 Fluorescence attachment, 6 position, UV shield; 100 W HBO mercury vapor lamp, with collimating lens, with external PSU BUT without objectives,

DX.9749 Empty filter cube for 6-position fluorescence attachment

without filter sets and empty cubes

DX.9930 3 W NeoLED™ replacement unit 100 W 24 Vac halogen replacement bulb 100 W HBO mercury vapor lamp

AE.5130 Universal Ø 23.2 mm tube adapter with built-in 2x lens for SLR photo camera with APS-C sensor. Needs T2 adapter

AE.5025 T2 ring for Nikon D SLR digital camera AE.5040 T2 ring for Canon EOS SLR digital camera

Microscope slides 76 x 26 mm, ground edges, 50 pieces Cover glasses 18 x 18 mm, thickness 0.13-0.17 mm, 100 pieces

Cover glasses 22 x 22 mm, thickness 0.13-0.17 mm, 100 pieces Lens cleaning paper, 100 sheets per pack

Immersion oil, n = 1.482 (25 ml)

PB.5274 Isopropyl alcohol 99% (200 ml)

PB.5275 Cleaning kit: lens cleaning fluid, lint free lens tissue, brush, air blower, cotton swabs



The Euromex Quality system is certified according to ISO 9001:2008 and supports our pursuit of continuous improvement and our on-going commitment to provide our world-wide customers

HEADQUARTERS

Euromex Microscopen bv

Papenkamp 20 6836 BD Arnhem The Netherlands Tel: +31 (0) 26 323 22 11 info@euromex.com

SALES EXPORT OFFICE

Euromex Microscopen Spain sl Carretera de Barcelona 88, Entresuelo Edificio Technomar 08302 Mataró, Spain Tel: +34 (0) 937 415 609 info@euromex.com